



INSTRUCTIONS

DP22/DP27

MICROSCOPE DIGITAL CAMERA

This instruction manual is for the Olympus microscope digital camera DP22/DP27.

To obtain optimum performance of this camera and to ensure the safety, study this manual thoroughly before operating the camera and keep it on hand during operation of the camera.

Keep this instruction manual in a safe place.

For details of products included in the configuration of this system, see page 16 of this instruction manual.

Optical Microscope Accessory

This product is applied with the requirements of standard IEC/EN61326-1 concerning electromagnetic compatibility.

- Immunity Applied to industrial and basic environment requirements.



In accordance with European Directive on Waste Electrical and Electronic Equipment, this symbol indicates that the product must not be disposed of as unsorted municipal waste, but should be collected separately.

Refer to your local Olympus distributor in EU for return and/or collection systems available in your country.

NOTE: This product has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the product is operated in a commercial environment. This product generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications.

Operation of this product in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

FCC WARNING: Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the product.

For Korea only

B급 기기 (가정용 방송통신기자재)

이 기기는 가정용(B급) 전자파적합기기로서 주로 가정에서 사용하는 것을 목적으로 하며, 모든 지역에서 사용할 수 있습니다.

Table of Contents

Introduction	1
Safety precautions	2
Handling precautions.....	5
1 System diagram.....	16
2 Nomenclature of respective portions	17
3 Assembly.....	18
1 Attaching camera head	18
2 Connecting interface cables	20
4 Installation of software.....	21
5 Outline of image acquisition procedures	22

6 Specifications.....	23
-----------------------	----

7 Troubleshooting.....	26
------------------------	----

■ Proper selection of the power supply cord.....	32
--	----

Introduction

With this microscope digital camera DP22/DP27 attached to the Olympus UIS2/UIS series microscope using the camera adapter, the image from the optical microscope can be acquired.

Using this microscope digital camera in combination with microscopes of other manufacturers may not obtain the complete optical performance.

Contents in this instruction manual

This Instruction Manual covers only the contents concerning the microscope digital camera DP22 and DP27. For operating procedures according to the connection status of the microscope digital camera, refer to the Instruction Manual according to the relevant connection status.

Safety precautions

If the product is used in a manner not specified by this manual, the safety of the user may be imperiled. In addition, the product may also be damaged. Always use the equipment according to this instruction manual.

The following symbols are used in this instruction manual.

 **CAUTION** : Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.

 **CAUTION** : Indicates a potentially hazardous situation which, if not avoided, may damage the equipment or other property.

© : Indicates commentary (for ease of operation and maintenance).

CAUTION - Installation of microscope -

Do not allow the height of the microscope higher than 1 m.

In order to prevent overturning, do not combine accessories to allow the height of the microscope higher than 1 m.

以上内容仅为本文档的试下载部分，为可阅读页数的一半内容。如要下载或阅读全文，请访问：<https://d.book118.com/058102112076006100>