

Instructions

DP75

Microscope digital camera

Note

This instruction manual is for DP75.

To ensure safety, and obtain optimum performance of this product, we recommend that you study this manual thoroughly before operating this product, and always keep this manual at hand when this product is in use.

Retain this instruction manual in an easily accessible place near the work desk for future reference.

For details of products included in the configuration of this system, see page 「System chart」 Page12.

Optical Microscope Accessory

A M 5 5 5 4

This product is applied with the requirements of standard IEC/EN61326-1 concerning electromagnetic compatibility.

- Immunity Industrial and basic electromagnetic environment

Emissions exceeding the level required by aforementioned standards may occur if this product is electrically connected to other equipment.

NOTE: This product has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the product is operated in a commercial environment. This product generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this product in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

FCC WARNING: Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the product.

FCC Supplier's Declaration of Conformity

Hereby declares that the product

Product name: Optical Microscope Accessory

Model Number: DP75

Conforms to the following specifications:

FCC Part 15, Subpart B, Section 15.107 and Section 15.109

Supplementary Information:

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

Responsible Party Name: Evident Scientific, Inc.

Address: 48 Woerd Ave Waltham, MA 02453, USA

Phone Number: 781-419-3900

In accordance with European Directive on Waste Electrical and Electronic Equipment, this symbol indicates that the product must not be disposed of as unsorted municipal waste, but should be collected separately. Refer to your local distributor in EU for return and/or collection systems available in your country.

For Korea only

Contents

1. Introduction	1
2. Safety Precautions	4
2.1 General	4
2.2 Installation	4
2.3 Electric shock prevention	4
2.4 Electrical safety	5
2.5 Injury prevention	5
2.6 Preventing burn	5
2.7 Safety-related labels on the product	6
2.8 Cautions and labels	6
2.9 Handling Instructions	6
2.10 Cyber security	6
3. Restrictions	7
3.1 Specifications subject to restrictions depending on the combination	7
3.1.1 Restriction condition	7
3.2 Specifications subject to restrictions depending on usage conditions	8
3.2.1 Cameras in general	8
3.2.2 Camera images	8
3.2.3 Image data	9
3.2.4 Position navigator	9
3.2.4.1 Prerequisite	9
3.2.4.2 Performance constraints	9
3.2.5 PC combination	11
4. System chart	12
5. Nomenclatures	13
6. Assembly	14
6.1 Attaching the IR cut filter insertion/removal knob	14
6.2 Mounting the camera head	15
6.3 Connecting the interface cables	17
6.4 Connecting external devices	17
6.5 Connecting AC adapter	18
7. Installing software	19
7.1 Software cellSens/PRECiV	19
7.2 Before installing software	19
7.3 The TWAIN driver DP2-TWAIN	19
7.4 Selecting Device	19

8. Summary of steps to acquire the images	21
9. IR cut filter insertion/removal function	23
10. External triggering function	24
11. Maintenance	25
11.1 Cleaning each part	25
11.2 Cleaning IR cut filter	26
12. Troubleshooting	27
13. Specifications	32
13.1 Basic specification table	32
13.2 Table of setting items by software	35
13.3 Operating environment	36
14. Proper selection of the power supply cord	37

1. Introduction

The Microscope Digital Camera DP75 is a product to acquire and process digital images, which is attached to our microscopes using a C mount camera adaptor made by our company.

The product will not work if a C mount adaptor made by other manufacturer or the product is attached to a microscope made by other company.

Cautions and confirmation before use

When you unpack the package, check all the included items referring to the list of accessories. If any components are missing or damaged, contact an Evident distributor.

Configuration of instruction manuals


Read all the instruction manuals supplied with the units you have purchased.

The following instruction manuals are produced for the units to be used with DP75.

Manual names	Main contents
DP75 (This instruction manual)	Request for safety, setup procedures, maintenance, troubleshooting, specifications, etc.
DP75 Quick Reference Guide for cellSens	Software specifications Operating procedures, functions, cautions, useful information, etc.
DP75 Quick Reference Guide for PRECiV	Software specifications Operating procedures, functions, cautions, useful information, etc.

How to read this instruction manual

This instruction manual is structured as follows.

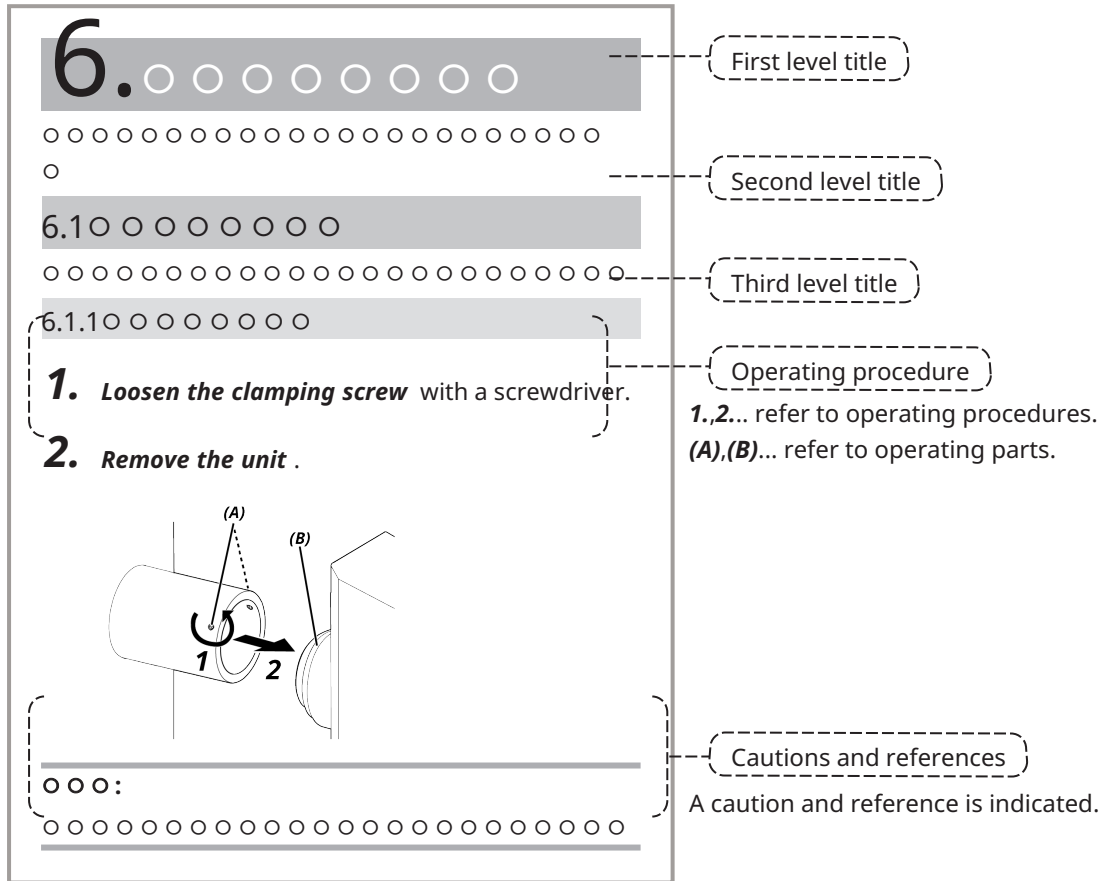

Cautions and references

This instruction manual is structured as follows.

⚠CAUTION:

Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury. It may also be used to alert the user to unsafe practices or potential material damage.

NOTE:

Indicates a potential situation which, if not avoided, may result in failure of this instrument.

TIP:

Indicates useful knowledge and information for use.

以上内容仅为本文档的试下载部分，为可阅读页数的一半内容。如要下载或阅读全文，请访问：<https://d.book118.com/466203242142010144>