



INSTRUCTIONS

DP26

MICROSCOPE DIGITAL CAMERA

This instruction manual is for the Olympus microscope digital camera DP26.
We recommend that you study this manual thoroughly before operating the camera and familiarize yourself fully with the use of this camera, which is important to ensure the safety, and to obtain optimum performance.

Retain this instruction manual in an easily accessible place near the work desk for future reference.

AX7970



This device complies with the requirements of directive 2004/108/EC concerning electromagnetic compatibility and of directive 2006/95/EC concerning electrical equipment designed for use within certain voltage limits. CE marking means the conformity to these directives.



In accordance with European Directive 2002/96/EC on Waste Electrical and Electronic Equipment, this symbol indicates that the product must not be disposed of as unsorted municipal waste, but should be collected separately.

Refer to your local Olympus distributor in EU for return and/or collection systems available in your country.

NOTE: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications.

Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

FCC WARNING: Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

For Korea only

B급 기기 (가정용 방송통신기자재)

이 기기는 가정용(B급) 전자파적합기기로서 주로 가정에서 사용하는 것을 목적으로 하며, 모든 지역에서 사용할 수 있습니다.

IMPORTANT – Be sure to read this section for safe use of the equipment. – 1-12

1	Intended use.....	4
2	Handling Precautions.....	5
3	Maintenance and Storage.....	6
4	Cautions for using.....	7
5	PC and Software	10

1	SYSTEM CHART	13
----------	---------------------------	-----------

2	NOMENCLATURE	15
----------	---------------------------	-----------

3 INSTALLATION..... 16

- 1** Installing the Camera Head 16
- 2** Connecting the Camera Cable 18
- 3** Connecting the AC adapter 19

4 IMAGE RECORDING PROCEDURE 21

- 4-1 Software 21
- 4-2 Image Recording Flow 22

5 SPECIFICATIONS 23

6 TROUBLESHOOTING GUIDE 25

IMPORTANT

The microscope digital camera DP26 is designed to be connected to a camera adapter mounted on an Olympus UIS2/UIS series microscope for recording of microscopic magnified images.

When the DP26 is used with a camera adapter from other manufacturers than Olympus, the optical performance may not be fully manifested.

Descriptions in Instruction Manual

This Instruction Manual contains only the descriptions of the microscope digital camera DP26. For procedures to use the connection kit of the microscope digital camera, refer to the Instruction Manual of the relevant connection kit.

If the equipment is used in a manner not specified by this manual, the safety of the user may be endangered. In addition, the equipment may also be damaged. Always use the equipment as outlined in this instruction manual.

The following symbols are used to set off text in this instruction manual.



: Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury or damage to the equipment or other property. It may also be used to alert against unsafe practices.



: Indicates commentary (for ease of operation and maintenance).

以上内容仅为本文档的试下载部分，为可阅读页数的一半内容。如要下载或阅读全文，请访问：<https://d.book118.com/666153124142010144>